**Bürkle GmbH**

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

## Safety instructions

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to escaping or off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood. Wear appropriate protective clothing.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ The instrument is suitable for indoor use and manual operation. Do not expose the device to any direct sunlight.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

### CAUTION!

Escaping liquids due to falling over or tipping of the instrument. Bodily injury due to escaping liquid and damage to the instrument.

- ▶ Make certain the container is positioned with adequate stability.
- ▶ Do not transport with the stopcock screwed on.

## Layout and function

- ▶ The device is made of PP, PE and PTFE.
- ▶ The device is suitable for dispensing non-crystallizing liquids (without solids) from canisters and barrels.
- ▶ The device can be used in the temperature range of 0 - 40 °C (briefly up to 60 °C).
- ▶ The device is pressure-resistant without a discharge up to max. 2 bar at 20 °C.

## Scope of delivery

1 x	Spigot	
1 x	Transparent LDPE seal	②

Point out any damage in the packaging to the freight forwarder/supplier immediately.

## Assembly and safe usage

### CAUTION!

Possible danger of breakage.

- ▶ Apply only normal manual force.
- ▶ Do not lift the canister by the stopcock.
- ▶ Do not suspend filling containers from the stopcock.

1. Insert the transparent LDPE seal ② in the sealing flange inside the union nut ①.
2. Screw the stop cock with union nut ① onto the container thread (G 3/4" connecting thread).
3. If applicable insert the discharge reducing nozzle ⑤.

### ⚠ WARNING!

Danger of bodily injury due to liquid under high pressure escaping.

- ▶ The pressure in the canister must not exceed 2 bar.
- ▶ When the discharge reducing nozzle ⑤ is being used, only a non-pressurized discharge is permitted (nozzle is only inserted).

### CAUTION!

Openings and spaces on the device. Danger of crushing.

- ▶ Be careful for spaces and openings when activating the plug.

4. Open the plug ⑥ (turn anti-clockwise) and dispense liquid.

### NOTE!

When filling, ensure sufficient adequate air ventilation of the container. If necessary, open the canister lid during the filling process or use a venting valve in screw cap, e.g. for Bürkle compact jerrycans (see accessories).

5. Close the plug (turn clockwise).
6. After filling, turn the angle discharge ④ upwards to prevent dripping. The union nut ③ must not be too tight for this, loosen the union nut ③ if necessary.

**⚠ WARNING!**

Bodily injury due to escaping liquid

- ▶ Close the stopcock completely after dispensing (set plug to at least „zero“).
- ▶ Keep the stopcock closed when it is not in use.
- ▶ Do not unscrew the stopcock from the container when it is filled.
- ▶ Empty the stopcock completely before disassembly.

**Troubleshooting**

Problem	To eliminate the error
Thread connection is not sealed	<ol style="list-style-type: none"> <li>1. Check transparent LDPE seal for position.</li> <li>2. Replace transparent LDPE seal with softer PE foam gasket (white) (see accessories).</li> <li>3. Thread is damaged (contact dealer or manufacturer).</li> </ol>
Stopcock is dripping	<ol style="list-style-type: none"> <li>1. Close the plug more tightly.</li> <li>2. Turn the discharge so it is pointing up.</li> </ol>

**Accessories**

Type	Specification	Number	Item No.
Compact jerrycan	5 l, with threaded connector 3/4"	1	0435-2050
Compact jerrycan	10 l, with threaded connector 3/4"	1	0435-1100
PE foam gasket	white	1	0006-5055
Venting valve in screw cap	PP/PE	1	0435-9200

More accessories can be found at [www.buerkle.de](http://www.buerkle.de)

**NOTE!**

When used with Bürkle compact jerrycan (see accessories), remove the transparent LDPE seal from the sealing flange of the union nut and insert the PE foam gasket (white) supplied with the compact jerrycan into the sealing flange inside the union nut.

- ▶ The compact jerrycan is suitable to fill off fluid and viscous liquids up to 500 mPas, pay attention to the chemical resistance.
- ▶ Constructive changes to the instrument or accessories are not permitted. Do not use PTFE thread sealing tape.

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