Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.

! WARNING!

The instrument is not suitable for highly flammable or explosive liquids.

- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Safety instructions during and after use

- ▶ The instrument generates overpressure. Possible bodily injury due to escaping liquids.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

Layout and function

- ▶ The disposable riser pipes are made of PE and the universal bungs of PVC.
- ▶ The instrument is suitable for filling thin to low viscosity liquids up to 500 mPas.
- ▶ The instrument is suitable at a temperature range-


from 0 to 40 °C.

! WARNING!

The system is not tight-closing. Possible gas leakage. Bodily injury due to escaping gases.

- ▶ Always keep valves open when not in use.
- ▶ Ensure adequate ventilation.
- ▶ Wear the necessary protective clothing.

- ▶ The bung fits onto container openings with a diameter of 30 to 70 mm forming an airtight seal.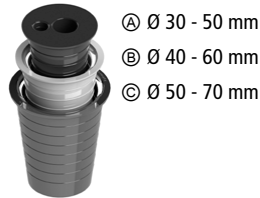

Ⓐ Ø 30 - 50 mm

Ⓑ Ø 40 - 60 mm

Ⓒ Ø 50 - 70 mm

Scope of delivery

10 x	Delivery tube, disposable	①
1 x	Universal bungs, three parts	②
1 x	Foot bellows	⑤
1 x	Cutter	③

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

1. Insert the delivery tube ① into the universal bung ② (opening "Rohr 12 mm") and push it through to the outlet.

NOTE!

If necessary adjust the universal bung to the opening diameter of the container.

- ▶ When using all three parts (Ⓐ, Ⓑ and Ⓒ) loosen Ⓑ and Ⓒ a bit of part Ⓐ to push the delivery tube through.

2. Determine the conveying tube length on the basis of the container height and shorten the delivery tube ① slightly at an angle to the container bottom using the cutter ③.

NOTE!

If necessary, shorten the discharge of the delivery tube ① with the cutter ③ to the desired length.


3. Insert the delivery tube ① with the universal bung ② into the container and press it in tightly.
4. Insert the stop-valve ④ into the universal bung ② (opening "Pumpe").

! WARNING!

Spraying of liquid under high pressure.

- ▶ Hoses and connections must be faultless.
- ▶ Ensure correct installation.

5. Close the stop-valve ④ (push the lever away from the connecting hose).

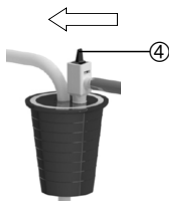**Safe filling**

1. Hold the discharge tube above the filling container.
2. Activate the foot bellows ⑤ and generate overpressure. Liquid starts to flow.
3. To complete the filling process open the stop-valve ④ (push the lever to the connecting hose) and allow the overpressure to escape.


Troubleshooting

Problem	To eliminate the error
The instrument does not generate an overpressure	<ol style="list-style-type: none"> 1. Stop-valve is opened. 2. The instrument is not connected with the container to form a seal.
Instrument is not pumping any liquid	<ol style="list-style-type: none"> 1. The instrument is not connected with the container to form a seal. 2. The riser pipe does not extend into the liquid being filled. 3. The delivery tube is not cut at an angle at the height of the tank bottom.

Accessories

Type	Specification	Number	Item No.
Delivery tube PE, disposable	For OTAL® disposable pumps	10	5000-8000

More accessories can be found at www.buerkle.de

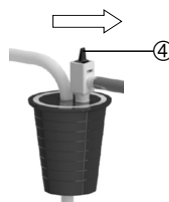**! WARNING!**

The container is under pressure. Bodily injury due to escaping liquid.

- ▶ Allow excess air to escape completely from the container.
- ▶ After pressure reduction, drain off any remaining liquid of the discharge tube completely.

NOTE!

The stop-valve is not suitable for precise dosing.

Completing filling and dismantling

1. Pull the stop-valve ④ off the universal bung ②.
2. Carefully remove the universal bung ② with delivery tube ① from the container.
3. Cut off the delivery tube ① directly below the universal bung ② using a cutter ③.
4. Pull the upper end of the delivery tube ① out of the universal bung ② and dispose it properly with the lower end.

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