Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ Do not fill the instrument with potentially explosive substances.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ Do not use any force when operating the instrument.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

Safety instructions during and after use

NOTE!

Abrasion of the cutting edge of the boring tube ②.

- ▶ Carry out a visual inspection every time before using.

- ▶ The instrument is suitable at a temperature range from 0 to 60 °C.
- ▶ Pay attention to have sufficient space above during the application (minimum height: double length of the sampler).

Layout and function

- ▶ The instrument is made of stainless steel V4A.

- ▶ The instrument is used to take samples of bulk materials which, due to their particle form, can easily become stuck or settle during sampling.
- ▶ The device is suitable for indoor and outdoor use.
- ▶ The insertion depth depends on the density of the bulk goods.
- ▶ The instrument is maintenance-free and suitable for steam sterilisation (bis 121°C).

Scope of delivery

1 x	Core sampler (consisting of boring rod ① and boring tube ②)	
1 x	Fixing clip	④

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

NOTE!

Bending of the boring rod ①.

- ▶ Do not plunge the instrument into highly compressed material.
- ▶ Do not perform any lever movements.

CAUTION!

Pointed parts on the instrument. Risk of injury.

- ▶ Do not pierce the instrument completely through the container.
- ▶ Keep persons and body parts out of the danger area.

CAUTION!

Entanglement of the sampler by moving parts in the machines.

- ▶ Do not use the instrument, if moving parts in the machine are operating.

Safe sampling

1. Insert the boring rod ① into the sample medium.
2. Put the boring tube ② over the boring rod ① and insert it into the sample medium by rotating it.
3. Insert the boring tube ② up to the centering ring ③ of the boring rod ①.
4. Hold the sampler at the boring rod ① and pull it out of the medium to be sampled.
5. Open the instrument and empty the sample on an appropriate surface or suitable collecting container.
6. Clean the instrument. Use the cleaning brush (see accessories) for the inside of tubes.

NOTE!

Optionally fix the boring tube ② with the fixing clip ④ fixieren (see picture 1), e.g. for horizontal sampling, for transport or storage.

NOTE!

Possible loss of boring rod ① during use. Possible loss of fixing clip ④ after use.

- ▶ After piercing the medium to be sampled, immediately put the boring tube ② over it and pull out the sampler on the boring rod ①.
- ▶ Keep the fixing clip ④ away from the medium to be sampled.

Accessories

Type	Specification	Number	Item No.
Cleaning brush	PVC/stainless steel, 40 cm	1	5304-0040
Cleaning brush	PVC/stainless steel, 80 cm	1	5304-0080
Sample beaker with cover (PP, sterilisable)	125 ml	1	7103-0125
Sampling bag SteriBag	60 ml	500	5344-0061
Sampling bag SteriBag StandUp	400 ml	500	5344-0400
Sampling bag SteriBag Blue	650 ml	500	5344-7008
SteriBag Premium tamper-proof sample bag	300 ml	100	5344-8000

More accessories can be found at www.buerkle.de

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