**Bürkle GmbH**

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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### General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

### Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ Do not fill the instrument with potentially explosive substances.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.

#### **WARNING!**

Damage to the respiratory tract and other organs due to escaping media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

#### **NOTE!**

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/disinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

### Layout and function

The instrument is used for removing pourable powders and granulates. It is made of stainless steel and polypropylene (PP). The provided sampling bottles are made of polyethylene (LDPE). The instrument is maintenance-free and suitable for steam sterilisation (up to 121°C).

### Parts in contact with medium

- ▶ Stainless steel 1.4404/AISI 316 L
- ▶ Polypropylene (PP)
- ▶ Polyethylene (LDPE)

### Scope of delivery

1 x	QuickPicker	①
2 x	Sampling bottle (LDPE, 250 ml)	②
1 x	Cleaning brush	
1 x	Directions for use	

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

### Safe sampling

#### **CAUTION!**

Openings with sharp edges on the instrument. Danger of cutting.

- ▶ Keep hands well away from openings with sharp edges.

1. Screw the sampling bottle ② onto the thread of the sampler ③.
2. Pierce the instrument into the powder/granulate up to the desired depth.

#### **WARNING!**

The container can be weakened and may break by plunging into it.

- ▶ Do not plunge the instrument into a pressurized container.

#### **NOTE!**

The instrument is designed for manual operation.

- ▶ Do not hammer the instrument into the container by using a hammer.

3. The instrument is filled with the sample.
4. The sample is emptied through the open end into the sampling bottle.
5. Remove the instrument from the powder/granulate.
6. Seal the puncture at the container.
7. Clean the instrument. Unscrew the tip and use the provided cleaning brushes for the inside of the tube.

**CAUTION!**

Pointed parts on the instrument. Risk of injury.

- ▶ Do not transport the instrument when the tip is not secured. Do not throw the instrument around.
- ▶ Do not pierce the instrument completely through the container.

**Accessories**

Type	Specification	Item No.
Sampling bottle	LDPE, 250 ml	0318-0250
Sampling bottle	PP, 250 ml	0319-0250
Transport case	for QuickPicker	3600-3000
Control seal close-it and close-it food & pharma		
More information and more accessories can be found at <a href="http://www.buerkle.de">www.buerkle.de</a>		

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