**Bürkle GmbH**

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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### General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

### Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument.

#### NOTE!

- Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.
- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
  - ▶ Pay attention to the material resistance.

### Safety instructions during and after use

- ▶ The instrument is not suitable for the use with wood, stone and metal.
- ▶ Do not use the instrument on live humans or animals.
- ▶ The instrument is not suitable for the use with stationary machine tools.
- ▶ Wear the appropriate protective clothing.

### Layout and function

- ▶ The instrument is made of stainless steel 1.4404 (AISI 316 L) / 1.4571 (AISI 316 Ti)
- ▶ The instrument is suitable to take samples from deep-frozen and semi-solid substances in the food and beverage industry by using hand-guided drills

with speed control.

- ▶ The instrument is maintenance-free and suitable for steam sterilisation (up to 121°C) for sterilization by flame treatment.

### Parts in contact with medium

- ▶ Stainless steel 1.4404 (AISI 316 L) / 1.4571 (AISI 316 Ti)

### Scope of delivery

1 x	Borer	①
1 x	Holder	②
1 x	Ejector	③
1 x	Directions for use	

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

### Assembly

1. Mount the holder ② on the drill.
2. Mount the borer ① onto the holder.

#### CAUTION!

- Moving parts on the instrument. Risk of crushing.
- ▶ Mount the holder carefully on the drill.
  - ▶ Mount the borer carefully onto the holder.

#### NOTE!

- Verify that the borer is attached correctly before use.

### Safe sampling

#### ⚠ WARNING!

- Bodily injury due to accidental contact with electrical power. Danger of electric shock.
- ▶ Keep the cable of the drill out of the cutting area.

#### ⚠ WARNING!

- Sharp parts on the instrument. Risk of injury.
- ▶ Keep hands well away from the cutting area.

**⚠ WARNING!**

Bodily injury due to drilling through the sampling material without adequate securing.

- ▶ Use a suitable base (e.g. cutting board).
- ▶ Use a suitable locking mechanism or secure the sampling material against twisting and shifting.

**⚠ WARNING!**

Bodily injury due to entanglement of body parts, hairs or clothes.

- ▶ Keep hands well away from the cutting area.
- ▶ Use a hairnet if necessary.
- ▶ Ensure to wear tight-fitting clothes.

**⚠ WARNING!**

Recoil of the drill due to blocking of the borer.

- ▶ Secure the drill against twisting.
- ▶ Ensure to have enough support and stability.

**CAUTION!**

Moving parts on the instrument. Risk of crushing.

- ▶ Keep hands well away from moving parts.
- ▶ Place the borer carefully onto the sampling material.

**CAUTION!**

Risk of slipping and falling due to wet or smooth floor.

- ▶ Ensure to have enough support and stability.

**CAUTION!**

Break out of the cutting edges during use.

- ▶ Ensure not to drill on other materials than the sampling material.

**CAUTION!**

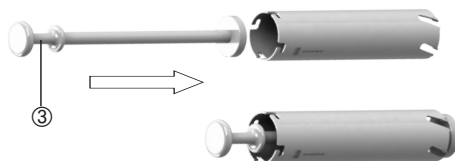
Frostbites due to contact with deep-frozen sampling material.

- ▶ Minimize contact time.

1. Place the borer onto the sampling material.
2. Start boring slowly until a slight trace is incised in the sampling material.
3. Accelerate the borer carefully.
4. Drill the borer to the desired depth into the sampling material.
5. Pull the borer out of the sampling material.

**After taking the sample**

1. Dismount the borer from the holder.
2. Empty the sample by using the ejector ③ into a suitable collecting container.

**CAUTION!**

Moving parts on the instrument. Risk of crushing when emptying the sample.

- ▶ Empty the sample carefully by using the ejector.

3. Dismount the holder from the drill.
4. Clean the instrument.

**CAUTION!**

Sharp parts on the instrument. Risk of injury.

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