Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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English

General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ Damage to the respiratory tract and other organs due to off-gassing media. Use in ventilated rooms or under a suitable exhaust hood.
- ▶ This instrument is not suitable for sampling highly flammable media. Not suitable for EX areas.

Safety instructions during and after use

- ▶ The instrument must only be used within the confines of applicable material resistance and for its

designated purpose. Any other type of use will affect the warranty/guarantee.

- ▶ Protect the device from moisture.
- ▶ Do not store the instrument with filled sample bottle.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

NOTE!

Follow the cleaning and maintenance instructions for the Bosch drive unit and pump head.

Layout and function

- ▶ The instrument is suitable for sampling liquids (max.

500 mPas) from a maximum suction depth of 5 m.

- ▶ The instrument is suitable for sampling liquids from 0 - 40 °C.
- ▶ For chemically demanding applications use pump hose made of Marprene (beige, durable), for food applications use pump hose made of silicone (transparent).
- ▶ The instrument is only suitable to take samples and not for continuous operation.
- ▶ The instrument is suitable for filling open sampling bottles.
- ▶ Only use the instrument with the recommended sample bottles specified in the accessories.
- ▶ Protect the instrument against effects of weather exposure during storage.

Parts in contact with medium

- ▶ PE (Suction hose)
- ▶ PP (Hose connector)
- ▶ Marprene or silicone (Pump hose)
- ▶ LDPE (Sample bottle)


Scope of delivery

2 x	Pump hose (1x Marprene, 30 cm; 1 x Silicone 30 cm)	①
1 x	Pump head	②
1 x	Suction hose , PE, 5 m	③
1 x	Hose connector, PP	④
1 x	Bottle basket	⑤
1 x	Sample bottle, LDPE, 250 ml	⑦
1 x	Bosch drive unit	⑩
2 x	Rechargeable battery	⑭
1 x	Battery charger	⑮

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

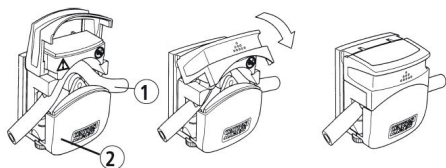1. Insert the pump hose ① into the pump head ②.

2. Cut the suction hose ③ off to the required length and attach it to the pump hose ① with the hose connector ④.
3. Fix the bottle basket ⑤ in place with the tightener ⑥.

CAUTION!

The mounting bracket may break.

- ▶ Do not load the suspension too heavily (max. 2 kg).
- ▶ Check the bottle basket to ensure it is properly seated. Allow the end of the holding rod to protrude out about 5 cm upward.
- ▶ Wear appropriate protective clothing.

4. Place the sample bottle ⑦ in the bottle basket ⑤.
5. Insert the pump hose ① about 1 - 3 cm into the sample bottle ⑦.

NOTE!

Allow the pump hose to protrude loosely into the sample bottle.

CAUTION!

Bursting of hoses and hose connectors due to wear or chemical stress.

- ▶ Carry out a visual inspection every time before using.

6. Place the pump attachment ⑧ in the holder ⑨ of the Bosch drive unit ⑩. Turn the locking ring ⑪ in the direction  until you hear it snap into place.

NOTE!

Follow the safety instructions for the Bosch drive unit and pump head.

Safe sampling

CAUTION!

Possibility of electrical shock.

- ▶ Follow the instructions of the drive manufacturer/battery charger manufacturer.

1. Set the gear selector switch ⑬ to „Gear 1 “ (low speed range).
2. Lower the suction hose ③ into the sample liquid.
3. Press the On/Off switch ⑫ to take the sample.

CAUTION!

Risk of atomisation. Ingress of liquids in the instrument.

- ▶ In case of uncontrolled escape of hazardous liquid, interrupt the sampling immediately and remove the Bosch drive unit.
- ▶ Contact the dealer or manufacturer.

NOTE!

For higher pumping heights it may take a while for the hose to be completely filled.

! WARNING!

Bodily injury due to escaping media if the sample bottle is overfilled.

- ▶ Do not overfill the sample bottle.
- ▶ Wear appropriate protective clothing.

4. To end sampling, release the On/Off button ④.

After taking the sample

NOTE!

Do not set the instrument down with the sample bottle filled. Remove the Bosch drive unit before setting it down.

1. Turn the pump head ② in the direction of  and pull it off the holder ③.
2. Remove the pump hose ① from the pump head ②.
3. Take the sample bottle ⑦ out of the bottle basket ⑤ and close it.

! WARNING!

Personal injury due to escaping media if the sample bottle is removed and the suction hose is taken out of the instrument.

- ▶ Disassemble the instrument carefully.
- ▶ Wear appropriate protective clothing.

4. Loosen the bottle basket ⑤ with the tightener ⑥.
5. If necessary rinse the hoses and hose connector with a pH-neutral liquid.

Accessories

Type	Specification	Number	Item No.
Suction hose Marprene for Vampire	100 cm	1	5327-0508
Suction hose Silicone for Vampire	5 m	1	8803-0609
Suction hose Silicone for Vampire	25 m	1	8881-0609
Sample bottle LDPE	250 ml	1	0318-0250
Sample bottle LDPE	500 ml	1	0318-0500
Sample bottle LDPE	1000 ml	1	0318-1000
Suction hose PE on roll 6 x 8 mm	Roll with 10 m	1	8805-0608
Suction hose PE on roll 6 x 8 mm	Roll with 100 m	1	8878-0608
Basket for sample bottle 250 ml	Ø 68 mm	1	5302-2025
Basket for sample bottle 500 ml	Ø 84 mm	1	5302-2051
Basket for sample bottle 1000 ml	Ø 100 mm	1	5302-2101

More accessories can be found at www.buerkle.de

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