Bürkle GmbH

Rheinauenstr. 5 | 79415 Bad Bellingen
Telefon +49 7635 6121-0
info@buerkle.de
www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) – and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument.

NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Pay attention to the material resistance.

Safety instructions during and after use

- ▶ The instrument is not suitable for the use with wood, stone and metal.
- ▶ The instrument is not suitable for the use with stationary machine tools.
- ▶ Wear the necessary protective clothing.
- ▶ The instrument is not suitable for cleaning in a dishwasher.

Layout and function

- ▶ The instrument is made of stainless steel 1.4034, 1.4301 and aluminium.
- ▶ The instrument is suitable for taking samples from all types of stalk material (e.g. hay, silage, straw, herbs, shrubs) using hand-held drills with speed control.
- ▶ The instrument is maintenance-free and suitable for steam sterilisation (up to 121 °C) or sterilisation by flaming.

Parts in contact with medium

- ▶ Drill: Stainless steel 1.4034 and aluminium
- ▶ Ejector: stainless steel 1.4301

Scope of delivery

1 x	Sampling drill collector tube	①
1 x	Sampling drill holder	②
1 x	Sampling drill drilling tip	③
1 x	Ejector	④

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

1. Mount the holder ② on the drilling machine.
2. Mount the collection tube ① on the holder ②.
3. Carefully mount the drill tip ③ on the collection tube.

CAUTION!

Moving parts on the instrument. Risk of crushing.

- ▶ Mount the holder carefully on the drilling machine.
- ▶ Carefully assemble the collection tube and drill tip.

NOTE!

Bodily injury due to ejected parts.
Check the fit of the drill before use.

⚠ WARNING!

Bodily injury due to accidental contact with electrical power. Danger of electric shock.

- ▶ Keep the cable of the drilling machine out of the cutting area.

⚠ WARNING!

Sharp parts on the instrument. Risk of injury.

- ▶ Keep hands well away from the cutting area.

Safe sampling

WARNING!

Bodily injury due to drilling through the sampling material without adequate securing.

- ▶ Use a suitable base (e.g. cutting board).
- ▶ Use a suitable locking mechanism or secure the sampling material against twisting and shifting.

WARNING!

Bodily injury due to entanglement of body parts, hairs or clothes.

- ▶ Use a hairnet if necessary.
- ▶ Ensure to wear tight-fitting clothes.

WARNING!

Recoil of the drilling machine due to blocking of the drill.

- ▶ Secure the drilling machine against twisting.
- ▶ Ensure to have enough support and stability.

CAUTION!

Moving parts on the instrument. Risk of crushing.

- ▶ Keep hands well away from moving parts.
- ▶ Place the borer carefully onto the sampling material.

CAUTION!

Risk of slipping and falling due to wet or smooth floor.

- ▶ Ensure to have enough support and stability.

CAUTION!

▶ Fire hazard! Heat development due to friction in the sampling material.

- ▶ Observe the temperature development in the sampling material and the surroundings.

CAUTION!

Break out of the cutting edges during use.

- ▶ Ensure not to drill on other materials than the sampling material.

1. Place the drill onto the sampling material.
2. Start drilling slowly. (Direction of rotation is clockwise)
3. Accelerate the drill carefully.
4. Drill into the sampling material to the desired depth.
5. Pull the drill out of the sampling material.

After sampling

1. Dismount the collection tube from the drilling machine.
2. Carefully dismantle the drill tip from the manifold.
3. Empty the sample by using the ejector into a suitable collecting container.
4. Dismount the holder from the drilling machine.
5. Clean the instrument carefully.

CAUTION!

Moving parts on the instrument. Risk of crushing when ejecting the sample.

- ▶ Empty the sample carefully by using the ejector.

CAUTION!

Sharp parts on the instrument. Risk of injury.

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