**Bürkle GmbH**

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

## Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ Do not fill the instrument with potentially explosive substances.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.

### **WARNING!**

Damage to the respiratory tract and other organs due to escaping media.

- ▶ When sampling hazardous substances, follow the safety data sheet of the medium to be sampled.
- ▶ Wear the appropriate protective clothing.

## LAYOUT and function

The instrument is used for removing pourable powders and granulates. It is made of stainless steel. The instrument is maintenance-free and suitable for steam sterilisation (121°C).

## Parts in contact with medium

- ▶ Stainless steel 1.4404/AISI 316 L

## Scope of delivery

|       |                                 |   |
|-------|---------------------------------|---|
| 1 x   | StickProof                      | ① |
| 1 x   | Clamp                           | ② |
| 100 x | PE sampling bags (120 x 170 mm) |   |
| 1 x   | Directions for use              |   |

- ▶ Check delivery for completeness and transport damage.
- ▶ Point out any damage in the packaging to the freight forwarder/supplier immediately.

## Safe sampling

### **CAUTION!**

Openings with sharp edges on the instrument. Danger of cutting.

- ▶ Keep hands well away from openings with sharp edges.

1. Attach the sampling bag with the clamp ② at the open end of the sampler ③.
2. Pierce the instrument into the powder/granulate up to the desired depth.

### **WARNING!**

The container can be weakened and may break by plunging into it.

- ▶ Do not plunge the instrument into a pressurized container.

### **NOTE!**

The instrument is designed for manual operation.

- ▶ Do not hammer the instrument into the container by using a hammer.

3. The instrument is filled with the sample.
4. The sample is emptied through the open end into the sampling bag.
5. Remove the instrument from the powder/granulate.
6. Seal the puncture at the container.
7. Clean the instrument.

### **CAUTION!**

Pointed parts on the instrument. Risk of injury.

- ▶ Do not transport the instrument when the tip is not secured. Do not throw the instrument around.
- ▶ Do not pierce the instrument completely through the container.

## Accessories

| Type   | Spezifikation        | Number | Item No.  |
|--|----------------------|--------|-----------|
| Clamp  | Stainless steel      | 1      | 5338-0001 |
| Packaging bags with closure  | 170 x 120 mm, 400 ml | 100    | 2348-0005 |
| Control seal close-it and close-it food & pharma   |                      |        |           |
| More information and more accessories can be found at <a href="http://www.buerkle.de">www.buerkle.de</a> |                      |        |           |

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