**Bürkle GmbH**

Rheinauen 5 | D - 79415 Bad Bellingen  
Telefon +49 (0) 76 35 8 27 95-0  
Telefax +49 (0) 76 35 8 27 95-31  
info@buerkle.de | www.buerkle.de

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## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.

## Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are sampling.
- ▶ When sampling hazardous substances, follow the technical rules on hazardous substances (in accordance with TRGS – (German) Technical Rules on Hazardous Substances) and wear appropriate protective equipment.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument.

### NOTE!

Before using, the instrument has to be cleaned to the extent of your (validated) cleaning process.

- ▶ Measures of manual cleaning have to be carried out in the cleaning/desinfection area and under consideration of occupational health and safety.
- ▶ Pay attention to the material resistance.

## Safety instructions during and after use

- ▶ The instrument can be used at temperatures up to 60 °C.
- ▶ The sampling area has to be free of disturbing factors and hazard sources (e.g. rotating elements).
- ▶ Do not perform any lever or batting movements during the application. The instrument may be deformed.
- ▶ Only operate the instrument with manual force.
- ▶ Pay attention to have sufficient space during use.
- ▶ The sampling area must be observable during the use.
- ▶ Wear the appropriate protective clothing.

## Layout and function

- ▶ The instrument is made of stainless steel 1.4404 (AISI 316 L).

- ▶ The instrument is suitable to take surface samples of frozen, highly viscous, paste-like food such as ice-cream and similar substances.
- ▶ The instrument is maintenance-free and suitable for steam sterilisation (up to 121°C).

## Parts in contact with medium

- ▶ Stainless steel 1.4404 (AISI 316 L)

Point out any damage in the packaging to the freight forwarder/supplier immediately.

## Safe sampling

### CAUTION!

Sharp parts on the instrument. Risk of injury.

- ▶ Keep hands well away from the cutting area.
- ▶ Do not insert the instrument into the mouth.

### CAUTION!

Frostbites due to contact with deep-frozen sampling material.

- ▶ Minimize contact time.
- ▶ Wear the appropriate protective clothing (cold protection gloves).

1. Connect the collection cylinder ① to the instrument using the bayonet connection ②.

### CAUTION!

Moving parts on the instrument. Risk of crushing.

- ▶ Attach the collection cylinder carefully.

### NOTE!

Possible loss of the collection cylinder during use.

- ▶ Verify that the collection cylinder is attached correctly before use.

2. Screw the instrument clockwise into the medium to be sampled.
3. Pull the instrument out of the medium to be sampled.
4. Disconnect the collection cylinder from the instrument.
5. Carefully remove the sample from the instrument.
6. Clean the instrument.

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