Bürkle GmbH

Rheinauen 5 | D - 79415 Bad Bellingen
Telefon +49 (0) 76 35 8 27 95-0
Telefax +49 (0) 76 35 8 27 95-31
info@buerkle.de | www.buerkle.de

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General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.
- ▶ The instrument is designed for use in industry, laboratories and science. The instrument is not suitable for any application in private or public spaces.

Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to escaping or off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood. Wear appropriate protective clothing.
- ▶ The instrument is not suitable for filling off combustible liquids.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument.

Safety instructions during and after use

CAUTION!

Escaping liquids due to falling over or tipping of the instrument. Bodily injury due to escaping liquid and damage to the instrument.

- ▶ Ensure sufficient stability of the instrument.
 - ▶ Shut down and secure the instrument immediately.
 - ▶ Contact the dealer or manufacturer.
- ▶ The instrument is suitable for indoor use and manual operation.
 - ▶ Pump with manual operation. Not suitable for continuous operation.
 - ▶ Pay attention to have sufficient space above during the filling process (minimum height: height of the maximum extended piston + 20 cm).
 - ▶ The instrument is not suitable for cleaning in a dishwasher.

NOTE!

Pump and valves can be damaged or plugged by dirt particles.

- ▶ Do not fill off any polluted media.
- ▶ Rinse the pump with a pH-neutral medium.

Layout and function

- ▶ The instrument is made of non-conductive material (PP, PVC) and is suitable for filling acids and alkalis (pay attention to the chemical resistance).
- ▶ The instrument is suitable for filling from containers with a capacity of 60 - 1000 l into smaller containers.
- ▶ The instrument can be used with a discharge tube with maximal length of 1,20 m.
- ▶ The instrument is maintenance-free and suitable at a temperature range from 0 to 40 °C.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

Assembly

⚠ WARNING!

Some liquid will continue to run while the discharge hose is being screwed on. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valve ③ on the discharge hose before mounting.
1. Screw the discharge hose ② onto the riser pipe ①.
 2. Guide the barrel screw connection (accessory; not included in delivery) and riser pipe ① with the discharge hose ② into the container and screw it together finger-tight.
 3. If necessary adjust the riser pipe.

NOTE!

Ensure to use a fitting thread (see accessories).

- ▶ Screw the thread by hand and without any force.

Safe filling

1. Hold the discharge hose in the filling container.
2. Open the shut-off valve ③ on the discharge hose.
3. Activate the piston rod ④ by using the handle ⑤ and fill off the liquid.

CAUTION!

Fluid ejection off the filling container. Bodily injury due to escaping liquid.

- ▶ Wear the necessary protective clothing.

⚠ WARNING!

Some liquid will continue to run after the discharge hose is unscrewed. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valve on the discharge hose.
- ▶ Raise the discharge hose and allow the liquid to run back into the container.

Completing filling and dismantling

1. After the last pump stroke, leave the piston rod in the lower position and insert it into the pump up to the stop.
2. Close the shut-off valve on the discharge hose.
3. To prevent any liquid continue to run (siphoning principle), fix the discharge hose above the liquid level in the original container.
4. Unscrew the discharge hose.
5. Carefully unscrew the instrument from the container.
6. Rinse the instrument with a pH-neutral fluid.

Accessories

Type	Specification	Number	Item No.
PP barrel screw joint	R 2", steel fine outer thread	1	5600-3130
PP barrel screw joint	Mauser 2", coarse outer thread	1	5600-3170
PP barrel screw joint	TriSure, coarse outer thread	1	5600-3180
Container screw joint	Opening Ø 60 mm	1	5600-3140
PVC barrel plug	Ø 40 - 70 mm	1	5600-3150
Discharge hose with stopcock	1,20 m	1	5600-0100
SafetyBox	Incl. safety goggles and protective gloves	1	2507-9000
Drain safety stand	Total length 117 cm, Ø 35 mm	1	5618-1000

More accessories can be found at www.buerkle.de

Troubleshooting

Problem	To eliminate the error
Liquid is not pumped.	Bottom valve is polluted. Clean the pump. If there is only little residual liquid in the container, the pumping must be carried out for a longer time.

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