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Directions for use | Instrucciones para el uso | Instruction d'emploi | Инструкция по пользованию

Art.-Nr. 5610-1030 / 5610-1040

### Fasspumpe Gasdicht PP

Gas-tight barrel pump PP

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## General safety requirements

- ▶ Read these directions for use carefully before using.
- ▶ Keep these directions for use where they will be accessible for all users.
- ▶ Failure to comply with the directions for use shall render all warranty conditions null and void and may lead to personal injury and material damage.
- ▶ The instrument must only be used by qualified skilled employees.
- ▶ The instrument is designed for use in industry, laboratories and science. The instrument is not suitable for any application in private or public spaces.

## Safety instructions before using

- ▶ Follow the manufacturer's instructions and safety instructions of the medium you are filling.
- ▶ Damage to the respiratory tract and other organs due to escaping or off-gassing media. Fill only in ventilated rooms or under a suitable exhaust hood. Wear appropriate protective clothing.
- ▶ The instrument is not suitable for filling off combustible liquids.
- ▶ The instrument must only be used within the confines of applicable material resistance and for its designated purpose. Any other type of use will affect the warranty/guarantee.
- ▶ Do not use any force when operating the instrument.

## Safety instructions during and after use

### CAUTION!

Escaping liquids due to falling over or tipping of the instrument. Bodily injury due to escaping liquid and damage to the instrument.

- ▶ Ensure sufficient stability of the instrument.
  - ▶ Shut down and secure the instrument immediately.
  - ▶ Contact the dealer or manufacturer.
- 
- ▶ The instrument is suitable for indoor use and manual operation.
  - ▶ Pump with manual operation. Not suitable for continuous operation.
  - ▶ Pay attention to have sufficient space above during the filling process (minimum height: height of the maximum extended piston + 20 cm).
  - ▶ The instrument is not suitable for cleaning in a dishwasher.

### NOTE!

Pump and valves can be damaged or plugged by dirt particles.

- ▶ Do not fill off any polluted media.
- ▶ Rinse the pump with a pH-neutral medium.

## Layout and function

- ▶ The instrument is made of non-conductive material (PP, PVC, HDPE) and is suitable for filling acids and alkalis (pay attention to the chemical resistance).
- ▶ The instrument is suitable for filling from containers with a capacity of 60 - 1000 l and a depth of 40 - 67 cm into smaller containers.
- ▶ The instrument is maintenance-free and suitable at a temperature range from 0 to 40 °C.

The barrel pump is available in two versions:

- ▶ 5610-1030: With universal adapter (30 - 70 mm) for a variable connection to the container.
- ▶ 5610-1040: With two thread adapters 3/4 „ with lock nut for a tightly screwed connection to the container.

Point out any damage in the packaging to the freight forwarder/supplier immediately.

## Assembly

### ⚠ WARNING!

Some liquid will continue to run. Bodily injury due to escaping liquid.

- ▶ Close the shut-off valves ③ on the discharge hose ① and the gas-shuttle pipe ② before mounting.

1. Pull out the telescopic riser pipe completely.
2. Press the bung of the pump ④ into the opening of the original container. Thereby, the telescopic riser pipe is adjusted to the ideal length.
3. Press the bung of the discharge hose ⑤ onto the filling container (version 5610-1030) or screw the thread adapters ⑥ onto the filling container (version 5610-1040).

### CAUTION!

Possible build up of dangerous overpressure.

- ▶ Drill holes in the bung must not be plugged.
- ▶ Do not kink the gas-shuttle pipe.

**NOTE!**

- Check the hose connections and seals regularly for position and tightness.

**Safe filling****CAUTION!**

Liquid entering the instrument. Bodily injury due to escaping liquid and damage to the instrument.

- Do not overfill the filling container.
- Do not fill off when the shut-off valves are closed.

1. Open the shut-off valves on the discharge hose and the gas-shuttle pipe.
2. Activate the piston rod ⑦ by using the handle ⑧ and fill off the liquid. Thereby, the displaced air is returned from the filling container to the original container via the gas-shuttle pipe.

**Completing filling and dismounting**

1. After the last pump stroke, leave the piston rod in the lower position and insert it into the pump up to the stop.
2. Carefully remove the bung from the filling container (version 5610-1030) or screw off the thread adapters (version 5610-1040) and allow the remaining liquid to drain out of the discharge hose into the filling container.
3. Close both shut-off valves.
4. To prevent any liquid continue to run (siphoning principle), fix the discharge hose above the liquid level in the original container.

** WARNING!**

Some liquid will continue to run. Bodily injury due to escaping liquid.

- Close the shut-off valves on the discharge tube and the gas-shuttle pipe before dismounting.
- Raise the discharge hose and allow the liquid to run back into the container.

5. Carefully remove the pump from the original container and allow the remaining liquid to drain into the filling container.
6. Allow the remaining vapours to escape out of the instrument and rinse the instrument with a pH-neutral liquid.

**Accessories**

Type	Specification	Number	Item No.
SafetyBox	Incl. safety goggles and protective gloves	1	2507-9000

More accessories can be found at [www.buerkle.de](http://www.buerkle.de)

**Troubleshooting**

Problem	To eliminate the error
Liquid is not pumped.	<p>Bottom valve is polluted. Clean the pump.</p> <p>The telescopic riser pipe does not extend into the liquid being filled.</p> <p>Open the shut-off valves.</p> <p>If there is only little residual liquid in the container, the pumping must be carried out for a longer time.</p>

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